

# Product manual

Labinco L31



LABINCO B.V.

P.O.BOX 3259-4800DG Breda, The Netherlands Zinkstraat 64-4823AC Breda, The Netherlands Tel.:0031(0)76-5417582, Fax:0031(0)76-5420173 E-mail:info@labinco-bv.com - Internet:www.labinco-bv.com



# Model L31



#### **Description:**

The Model "L31" is a rectangular hotplate 460x350x125mm, provided with a 2000 watt topplate for fast heat transfer.

Energy input control is provided, with stepless adjustable simmerstat controlled hotplate. Temperature from ambient up to 350 degr.C Cast steel coated housing.

## Operating instructions:

Install the unit on a flat sturdy bench in view of its weight and plug the power cord into a suitable current outlet. Take care that the heating plate does not stand too near to heat sensitive objects or the front of the bench. The heat regulator is very sensitive and it is therefore recommended, that once the best setting for a specific temperature and product load has been found that the unit be turned on or off with its illuminated power-switch. Because of the large power consumption it is recommended that the unit is put on a separate line and not together with other instruments on the same power outlet.

Specifications: 230V 2000watt 50/60HZ

<u>Dimensions:</u> 460x350x125mm weight; 11 KG



### LABINCO B.V.

P.O.BOX 3259-4800DGBreda,The Netherlands Zinkstraat 64-4823AC Breda,The Netherlands Tel.:0031(0)76-5417582,Fax:0031(0)76-5420173 E-mail:info@labinco-bv.com - Internet:www.labinco-bv.com